

# QuickTask™ Mop

## Flat mop with polyester/cellulose face

The QuickTask™ range consists of a variety of flat mops that can be easily attached to a choice of flat mop frames and handle assemblies.

QuickTask TASK0550 mop heads are constructed from a hydroentagled polyester/cellulose nonwoven fabric with a synthetic fiberfill core. This provides a highly sorbent mop which is suitable for the removal of light soil, residues and biofilms. Designed for single use, the mops are compatible with a wide range of chemicals and disinfectants.



Available irradiated.

| Features                                                  | Benefits                                                                                                                                                                                                                       |
|-----------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Highly sorbent fabric construction                        | <ul style="list-style-type: none"><li>• Allows for removal of light soils, residues, and biofilms</li><li>• Very absorbent and drip-free</li><li>• Compatible with most common solutions, solvents and disinfectants</li></ul> |
| Polyester/cellulose cover over a polyester fiberfill core | <ul style="list-style-type: none"><li>• Makes it ideal for cost-effective single use</li></ul>                                                                                                                                 |
| Available gamma irradiated                                | <ul style="list-style-type: none"><li>• Suitable for all grade of pharmaceutical cleanrooms</li></ul>                                                                                                                          |

| Part No. | Description                                         | Size                          | Packaging           |
|----------|-----------------------------------------------------|-------------------------------|---------------------|
| TASK0550 | QuickTask Mop<br>Irradiated, polyester/cellulose    | 16" x 5"<br>(406 x 130 mm)    | 1/bag; 40 bags/case |
| 2740     | QuickTask Mop Frame<br>Polypropylene                | 16" x 4.25"<br>(41 x 11 cm)   | 1/case              |
| 2699P    | QuickTask Magnetic Mop Frame<br>Polypropylene       | 15.5" x 3.75"<br>(40 x 10 cm) | 1/case              |
| 2906     | QuickTask Magnetic Mop Frame<br>Stainless steel     | 15.5" x 3.75"<br>(40 x 10 cm) | 1/case              |
| 6208     | QuickTask Non-magnetic Mop Frame<br>Stainless steel | 15.3" x 2.3"<br>(40 x 6 cm)   | 1/case              |

| Product Information   |                                                                                                                                                                                                                                                                                |
|-----------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Mop material          | • Nonwoven polyester/cellulose cover; Synthetic fiberfill core                                                                                                                                                                                                                 |
| Mop construction      | • Multilayer sewn lamination/sealed border                                                                                                                                                                                                                                     |
| Packaging materials   | • Outer bags (OB1, OB2, OB3), low density polyethylene (LDPE) <br>Case (CS), corrugated fiberboard (PAP)  |
| Environment           | • ISO 5-8    Grade A/B for sterile, C/D for nonsterile                                                                                                                                                                                                                         |
| Shelf life            | • Sterile: 2 years from manufacturing date                                                                                                                                                                                                                                     |
| Dimensions            |                                                                                                                                                                                                                                                                                |
| As packaged; nominal: | • 15.5" x 5.25" (39.4 x 13.3 cm)                                                                                                                                                                                                                                               |
| Mounted on frame:     | • 16" x 4.5" (41 x 11.5 cm)                                                                                                                                                                                                                                                    |



TASK0550

| Recycle Symbols                                                                     |                                                                                     |                                                                                     |                                                                                     |                                                                                      |
|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------|
| PET                                                                                 | HDPE                                                                                | LDPE                                                                                | PP                                                                                  | PAP                                                                                  |
|  |  |  |  |  |

| Technical Data                                  |                         |             |
|-------------------------------------------------|-------------------------|-------------|
| Attribute (units)                               | Typical Value           | Test Method |
| Mop weight, dry; (g)                            | 31                      | CTM004-03   |
| Sorbent capacity; (mL/mop)                      | 352                     | CTM004-03   |
| Particles ≥ 0.5µm; (x10 <sup>6</sup> /mop)      | 8.8                     | CTM006-01   |
| Fibers ≥ 100µm; (x 10 <sup>3</sup> /mop)        | 0.46                    | CTM006-01   |
| Mop coverage; ft <sup>2</sup> (m <sup>2</sup> ) |                         | CTM037-01   |
| S-Curve                                         | 347 (32)                |             |
| Place and Pull                                  | 224 (21)                |             |
| Mop frames                                      | 2906, 2699P, 2740, 6208 |             |

|          | EA/OB1 | OB1/OB2 | OB2/OB3 | OB3/CS | EA/CS |
|----------|--------|---------|---------|--------|-------|
| TASK0550 | 1      | 1       | 4       | 10     | 40    |

EA = each; OB = outer bag 1/2/3; CS = case

Notes

- a) The data shown are typical values and should not be used as product specifications.
- b) Valid product comparisons may only be obtained through side-by-side testing in the same test facility, under similar conditions.
- c) Current and/or comparison data may be available. Please contact a Contec sales representative for details.

Test Methods

- 1. CTM = Refers to Contec Test Method.
- 2. Mop coverage data is based on mop head applying solution such that the mopped surface is visibly wetted across the full mop path with no dry streaking observed, under controlled conditions. It is the responsibility of the user to determine appropriate coverage to ensure surfaces are visibly wet for the contact time required for their specific solution.